

cQrex[®] KC

Our next-generation cystine peptide to optimize biopharma cell culture

cQrex[®]



Chemically defined peptides for intensified bioprocesses


Efficient cell culture processes are the cornerstone of biological drug production. Evonik's cQrex® portfolio of chemically defined and highly pure peptides helps to formulate more concentrated and stable media, simplify the cell culture process, and increase productivity.

cQrex® KC is a highly soluble peptide designed to address the challenge of L-cystine supply in cell culture. It allows the formulation of concentrated and pH-neutral basal, feed and perfusion media, therefore intensifying and simplifying complex production processes. By increasing the cell proliferation, viability and cell-specific productivity, cQrex® KC also contributes to higher antibody titers and increased overall productivity of biological drugs.

Optimize your cell culture media with high concentrations of L-cystine

cQrex® KC is a precursor of the key amino acids L-cystine and L-lysine. The latter boosts solubility and is preferred over an alanine-based molecule when working with cells that are sensitive to L-alanine accumulation.

cQrex® KC (N,N'-di-L-Lysyl-L-Cystine dihydrochloride)

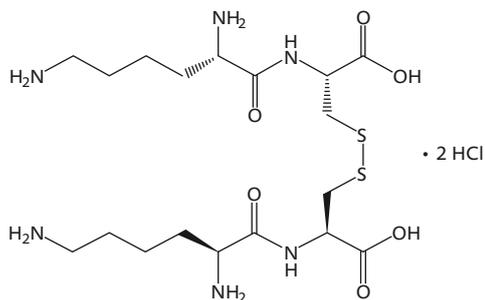

- Highly pure and chemically defined
- Efficiently metabolized by cells
- Non-animal derived
- Manufactured in-house in Europe


Want to discover more?

Get in touch with our experts:
cQrex@evonik.com

Overcome solubility limitations with cQrex® KC

- Solubility is around **1000 times higher than L-cystine** in pH-neutral media
- Rapid dissolution
- Increased flexibility in media formulation and bioprocess design

Max. solubility at neutral pH in water

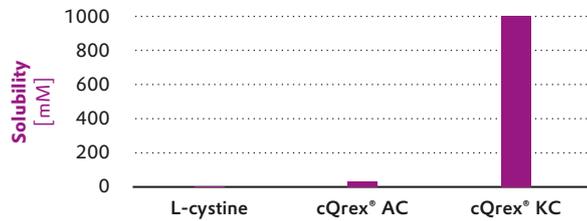


Maximize your bioreactor productivity with cQrex® KC

Fed-batch cultivation of a CHO-GS cell line expressing a monoclonal antibody in bioreactors:


- Equimolar amounts of L-cystine and cQrex® KC
- Adjusted L-lysine concentration

cQrex® KC supplementation in the feed has a positive impact on cell proliferation: it leads to a **prolonged high cell viability**.

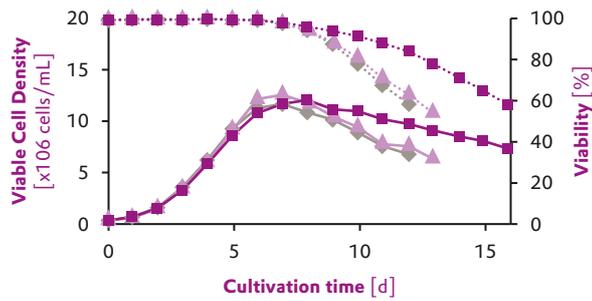

■ L-cystine in feed ■ cQrex® KC in feed ▲ cQrex® KC in basal medium

The longer cultivation time combined with a high cell-specific productivity enables a **strong increase in the antibody titer** when cQrex® KC is added to the feed or to the basal medium.

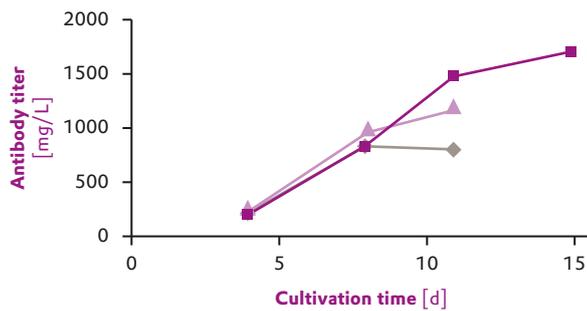

Cell specific productivity at d11

q_p [pg/(c*d)]

■ L-cystine in feed	10.4
■ cQrex® KC in feed	18.2
▲ cQrex® KC in basal medium	14.6

This information and all further technical advice are based on our present knowledge and experience. However, it implies no liability or other legal responsibility on our part, including with regard to existing third party intellectual property rights, especially patent rights. In particular, no warranty, whether express or implied, or guarantee of product properties in the legal sense is intended or implied. We reserve the right to make any changes according to technological progress or further developments. The customer is not released from the obligation to conduct careful inspection and testing of incoming goods. Performance of the product described herein should be verified by testing, which should be carried out only by qualified experts in the sole responsibility of a customer. Reference to trade names used by other companies is neither a recommendation, nor does it imply that similar products could not be used.

cQrex® – trademark of
Evonik Industries AG and its subsidiaries.

202302